

# Respiration Monitor Belt

The operation of the Respiration Monitor Belt requires that a **djb microtech** Low Pressure Sensor is connected to the belt. To make this easy, you have been supplied with two short lengths of Tygon tubing and two small white connectors.

To inflate your Respiration Monitor Belt carry out the following steps:

- Take the length of tubing that has the pointed connector and push the end of the tubing over the black pressure transducer that protrudes from the right hand side of the Low Pressure Sensor.
- Now connect the other length of tubing to the connector on the Respiration Monitor Belt.
- Push the two white connectors together.
- Turn the silver screw (near the hand pump) clockwise
- Now inflate the Respiration Monitor Belt.

Please note that the pressure sensor must be connected before inflating the belt.

## **djb microtech ltd**

Delfie House, 1 Delfie Drive, Greenock, Scotland, PA16 9EN

Phone 01475 786540 Email: [info@djb.co.uk](mailto:info@djb.co.uk) Website: [www.djb.co.uk](http://www.djb.co.uk)